

Brabender® Sedimat

Laboratory mill for the Zeleny
sedimentation test



ICC-Standard no. 116
ICC-Standard no. 118
ISO 5529

Brabender®

... where quality is measured.



Brabender Sedimat

For automatic production of test flours for the Zeleny sedimentation test.

The Sedimat mill, in combination with a shaker, is the instrumental prerequisite for the sedimentation test acc. to Zeleny (ICC Standard 116 and 118, ISO 5529).

This test describes the amount and quality of the gluten fraction of a flour. The sedimentation volume, determined in ml, provides information on the expected dough quality, gas retention capacity, fermentation tolerance and volume yield of the baked goods.

Yield, ash content and fineness of the test flours produced with the Sedimat meet the requirements for reproducible test performance.



Technical Data

Sedimat

- Capacity: approx. 100 g of wheat within 3 min.
- Mains connection:
3x 230 V; 50/60 Hz + PE; 2.0 A
3x 400 V; 50/60 Hz + N + PE; 1.1 A
- Dimensions (W x H x D): 380 x 620 x 610 mm
- Weight: approx. 50 kg net

Shaker

- Mains connection:
1x 230 V; 50/60 Hz + N + PE; 0.2 A
- Dimensions (W x H x D): 660 x 370 x 430 mm
- Weight: approx. 10 kg net



Highlights

Rolls made of high-quality special steel:

Prevent impairment of grinding results due to roll wear during normal operation.

Round sifter integrated in closed flour collecting vessel:


Excludes contamination of the flour sample to be milled with residues of the previous flour.

Shaker (optionally available):

Ensures homogeneous mixing with the chemicals and precise adherence to the proving and sedimentation times in accordance with Zeleny regulations.

Additional equipment for the Zeleny test:

- Shaker with up to 8 shaking cylinders
- Automatic dosing device

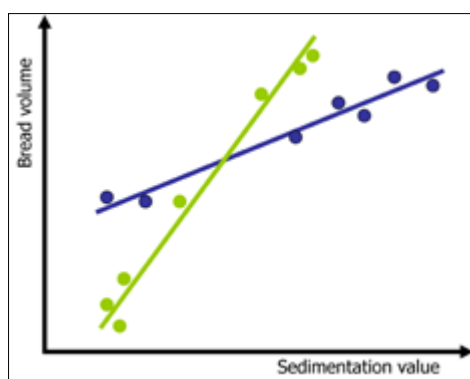

A Wheat: $420 + 3 \times \text{Sedi} + 10 \times \text{Protein} = \text{Bread volume}$
 B Wheat: $306 + 3 \times \text{Sedi} + 17 \times \text{Protein} = \text{Bread volume}$



Brabender® GmbH & Co. KG

Headquarter Germany
Kulturstraße 49 – 51
D-47055 Duisburg
Phone: +49 203 7788 0
sales@brabender.com
www.brabender.com

C.W. Brabender® Instruments, Inc.

USA
50 East Wesley Street
South Hackensack, NJ 07606
Phone: +1 201 343 8425
foodsales@cwbrabender.com
www.cwbrabender.com

000 Brabender®

Russia
Yagodinskaya St. 25
420032 Kazan
Phone: +7 843 233 46 66
000.brabender@brabender.ru
www.brabender.ru



Headquarter Germany



Contact us 24/7: +49 203 7788-131

Brabender agencies worldwide.
©2021 Brabender® GmbH & Co. KG

All trademarks are registered. Subject to design
and technical modifications without notice.

Brabender®
... where quality is measured.